


How to use our Magnetic Rack




Magnetic Separation Rack Instructions

Read before use

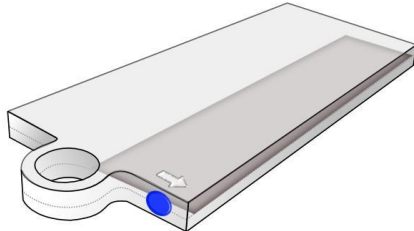


Magnetic insert


The colored dot and the arrow on the magnetic insert indicates the location of the magnetic field.

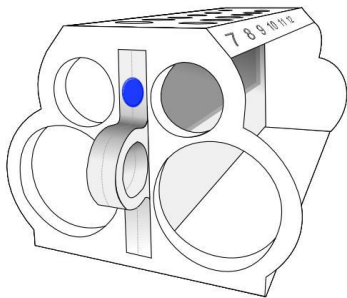

POSITION OF THE MAGNETIC FIELD
FOR 1.5mL or 15mL* TUBES

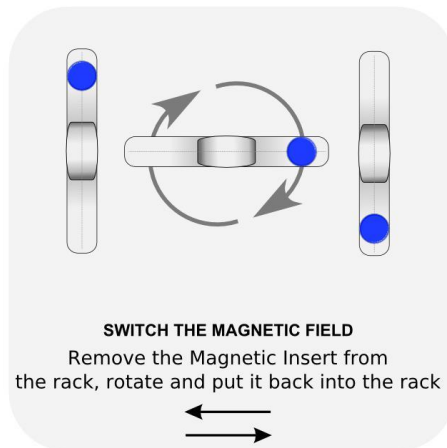

SWITCH THE MAGNETIC FIELD
Remove the Magnetic Insert from
the rack, rotate and put it back into the rack

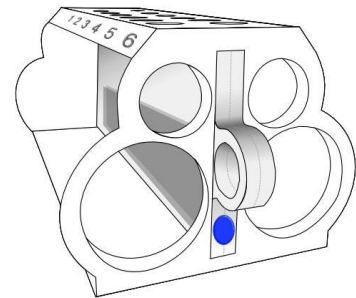

POSITION OF THE MAGNETIC FIELD
FOR 50mL* TUBES

* use it in vertical position.

Our Magnetic rack should be stored in a cool, dry environment and should be cleaned with a damp cloth and mild detergent or 70% ethanol. Do not autoclave. This magnetic rack should be handled with care, avoid direct contact with magnet material; pacemakers, implants, watches, electronic devices, all ferrous material and magnetic media have to be kept away.

We wish you the best for your experiments with OZ Biosciences' reagent.

OZ Biosciences' team

OZ Biosciences SAS
163 avenue de Luminy
Case 922, zone entreprise
13288 Marseille cedex 09 - FRANCE
Ph: +33 (0) 486 948 516
Fax: +33 (0) 463 740 015
contact@ozbiosciences.com
order@ozbiosciences.com

OZ Biosciences INC
7975 Dunbrook Road,
Suite B
San Diego CA 92126 - USA
Ph: + 1-858-246-7840
Fax: + 1-855-631-0626
contactUSA@ozbiosciences.com
orderUSA@ozbiosciences.com